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※For attendance confirmation, please write your full name and enrollment number on paper.

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Current Status of the Thin-section labo

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Thin-section room cleaning is carried out by users.

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薄片室 Thin-section labo

薄片室のルール Rules for Using the Thin-section Room

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The rules for using the thin-section room

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Available Hours:

Weekdays: 8:00 AM to 8:00 PM

Saturdays, Sundays, and Public Holidays: Not available for use

Records:

- When using the thin-section room, individuals must sign in and out in the entry management log (located at the door of the thin-section room)
- Record usage of rock cutters in the usage log.
- After use, cleaning is required
- For sound insulation purposes, entry and exit should be through the door of the precision polishing room.

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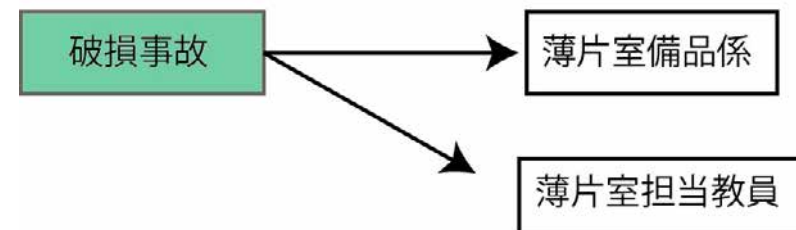
- ・ In case of an accident, immediately contact the designated personnel and prohibit the use of the affected equipment. Place a sign indicating the prohibition of use on the equipment in question.

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Injuries will be handled differently depending on their severity. In cases of severe injury, administer first aid such as hemostasis and immediately call for an ambulance (Tel: 119) to receive treatment at a medical facility.

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For minor cuts or injuries, such as a slight cut on the fingertip, use the first aid kit provided in the thin-section room for treatment.

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Before using any equipment, ensure there are no abnormalities:

Check all equipment for any signs of abnormalities (damage, unusual sounds, etc.).

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If any abnormalities are detected (such as damage or unusual noises), immediately contact the thin-section room supervisor (Professor Ueda).

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If abnormalities are found, do not use the equipment and place a notice prohibiting its use.

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**The rules for using machines in the
thin-section room**

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The rules for using machines in the thin-section room

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Always stabilize the specimen **using clay** before cutting.

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Inject grease into the bearing parts once a month to prevent wear and tear.

Fill any gaps with clay and firmly secure with both hands

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Attach the hose from the rear water tap and direct cooling water onto the blade.

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Insert the sample table placed aside into the cart.

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Secure the sample completely using clay.

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Turn on the main switch and operate by stepping on the foot switch.

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If the stone tilts during cutting and the blade locks, immediately turn off the power using the foot switch.

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After turning off the power, gently remove the stone and check if the blade is bent or chipped.

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Close the valve next to the blade to stop the water flow. Then, rinse away the rock debris with a sprinkler hose.

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Use the compressor to remove moisture from the stage.

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Wipe off any dirt that has splattered onto the walls or floor.

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Dispose of the rock debris on the desk and wipe it clean with a cloth.

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Before cleaning with water, turn off the power and elevate the foot switch above the floor to prevent electric shock.

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For details, refer to the notice posted next to the machine.

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After removing large rock debris by hand, allow mud and smaller debris to fall from the drain into the sediment trap (bucket) below.

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After use, wipe off any rock debris on the rails of the base with a cloth, then apply machine oil. Move the trolley back and forth to evenly distribute the oil.

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Lower the lid and lightly secure it with screws before use.

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It is acceptable to leave the blade attached.

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After use, rinse away any mud or debris inside using a sprinkler hose or sprayer.

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After use, rotate the blade and then wipe away moisture with a cloth.

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After use, apply KURE 556 to the slide rails.

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The particle size of the polishing powder used with the polishing disk (grinder) can be finer (higher number) than the specified size. (For example, it's acceptable to use 600 grit powder on a 120 grit polishing disk.)

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Black or green carborundum should be used in a dissolved state in water.

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If the polishing powder runs out, refill it from the bucket at your feet.

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After use, wipe the walls and tables with a cloth.

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Clean the grinder after use by rinsing off dirt with water and rotating it or blowing away moisture with an air duster.

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If water accumulates in the water receiver, pour off the excess into the sink.

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Widely used for tasks such as drying chips and cleaning equipment.

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When using an air duster, be cautious as the force can be strong and may cause chips or slide glasses to fly away.

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When using the sprayer, refill the sink's plastic bottle with water.

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During the weekend cleaning, turn off the main power, open the knob under the tank, and drain the accumulated water along with the air.

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Bringing and taking away chips during work is the basic practice.

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Dry and bond on aluminum foil with the student's grade, name, and time placed on it.

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Items that have been installed for more than 12 hours may be moved by individuals to the designated area for left chips, along with the aluminum foil.

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If there are no samples placed or only abandoned samples, the power should be turned off.

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Creating personal or laboratory-specific petropoxy from the shared one is prohibited.

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Petropoxy purchased by each laboratory should be stored within that laboratory.

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Petroproxy (clear) is typically mixed at a ratio of 10 (main agent) to 1 (hardener).

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Completely saturated silica gel (red) should be heated on a hot plate set at 120 degrees or in a constant temperature chamber until it turns blue.

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Cleaning duties are assigned to the person who used the equipment during that week.

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Every Friday from 18:00 onwards, the tasks are as follows:

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- ①Cleaning the hot plate
 - ②Draining the compressor
 - ③Cleaning the inside of the grinder and the water receiver
 - ④Cleaning the sediment trap
 - ⑤Floor cleaning
 - ⑥Disposing of garbage
 - ⑦Checking and replenishing consumables
- Tasks ③ and ④ are performed only at the end of the month.

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Every Friday Cleaning Tasks (For the Person Who Used the Equipment That Week):

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①Cleaning the Hot Plate:

- ・ Remove any left chips (it is acceptable to leave them overnight).
- ・ Remove any debris from inside the hot plate.
- ・ Use a spatula to scrape off hardened petro, smooth out any unevenness, and wipe with a wet cloth.

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②Draining the Compressor:

- ・ Turn off the compressor switch (press down the red button).
- ・ Place an aluminum tray under the tank, then twist and open the drain valve to release air and water.
- ・ Water may accumulate on the floor under the tank as it drains, so wipe it up with a cloth if necessary.
- ・ Once the sound of air being discharged stops and the pressure gauge indicating the air pressure reaches 0, close the valve and turn on the switch (pull up the red button).

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③Cleaning the Interior of the Polishing Disk and Water Receiver:

- ・ Follow the cleaning instructions outlined in the "Monthly Cleaning Tasks" section for the polishing disk.
- ・ Remove mud with a shovel.
- ・ Remove the urethane mat wrapped around the perimeter of the water receiver and wash it.
- ・ Dispose of the mud collected under the disk and the mud removed in step 2 at the designated stone disposal area and wash it.

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End-of-Month Cleaning Tasks (For the Person Who Used the Equipment That Month):

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④Cleaning the Sediment Trap:

Dispose of the accumulated mud in the sediment traps (white buckets) located under the sink, under the JAM cutter, and under the lower blade primary cutter at the designated stone disposal area. Wash the buckets with water from the stone disposal area's faucet.

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- Cleaning of Used Equipment
 - Rinse away any accumulated polishing agent in the sink.
 - Wipe down surfaces such as desks thoroughly.
 - Wash and then dry any used cloths.
- Let's keep everything clean for both the equipment and the next user!

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Combustible Waste:

- Kimwipes or newspaper
- Paper with dirt such as rock powder
- Cloth with dirt such as rock powder
- Paper with adhesive
- Paper used to wipe off lubricating oil
- Tapes
- Latex gloves
- Ethanol and acetone are volatile, so they can be treated as combustible waste.
- Remove moisture as much as possible.

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Plastic Waste:

- All plastic products
- Containers
- Plastic bags
- Teflon sheet scraps
- Styrofoam
- Saniment gloves
- Clumps of resin
- Syringes for chemical mixing

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Non-combustible Waste:

- Slide glass
- Slide glass with rock adhesive
- Polishing chips used with Specifics
- Waste containing a mixture of resin and stone
- Remove labels from metal cans and dispose of them as non-combustible waste.
- Aluminum foil
- Other metals

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※ Do not dispose of glass in the stone disposal area, as it may lead to injuries. This includes glass slides and glass that has been adhered to rocks.

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- Containers for chemicals → Waste liquid center
- Petro containers or rods
- Wipe syringes for reuse. If they are no longer usable, dispose of them as plastic waste.
- Remove labels from reagent bottles → Waste liquid center

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Items in the thin-section room are for communal use, so taking them out or keeping them is prohibited. If there is a valid reason for taking something out, consult the administrators or instructors.

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Necessary or commonly used items may be purchasable within the budget, so consider consulting to see if they can be acquired.

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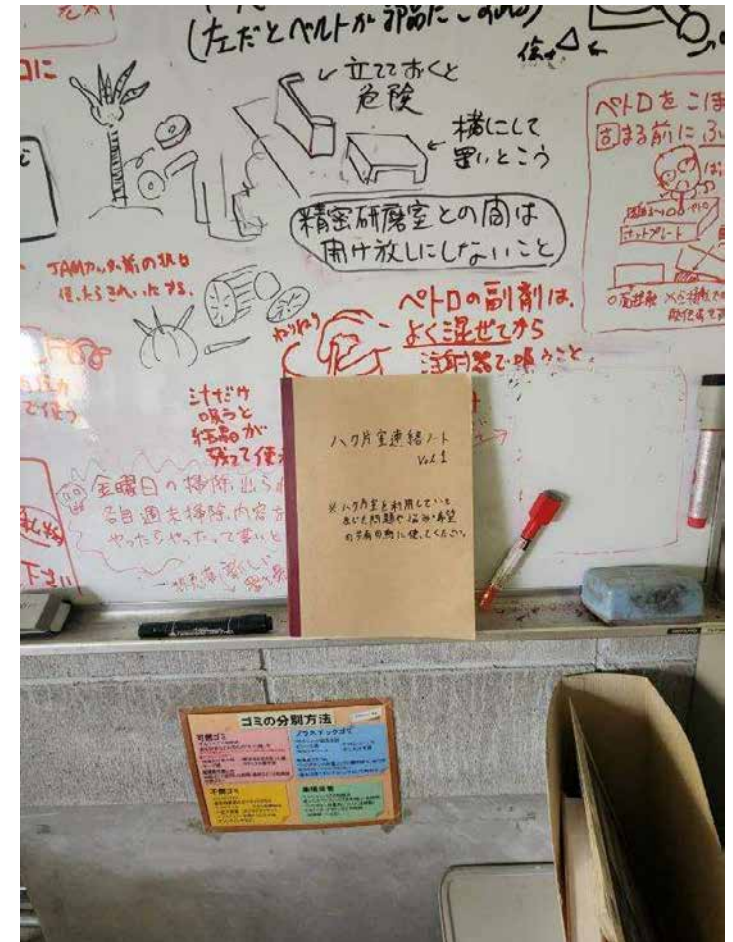
The thin-section room is a laboratory, and the equipment installed there is essential for everyone's research. Let's not forget to express our gratitude to the tools and equipment that work for us by thoroughly cleaning them after use and avoiding rough handling.

If you have any questions about the thin-section room, please feel free to ask the thin-section staff or Mr. Ueda.